Affyringsmekanismen består af

- Spole 1600 vindinger (462530)
- Jernkerne med fingerskrue
- Anker med stikben (195590)

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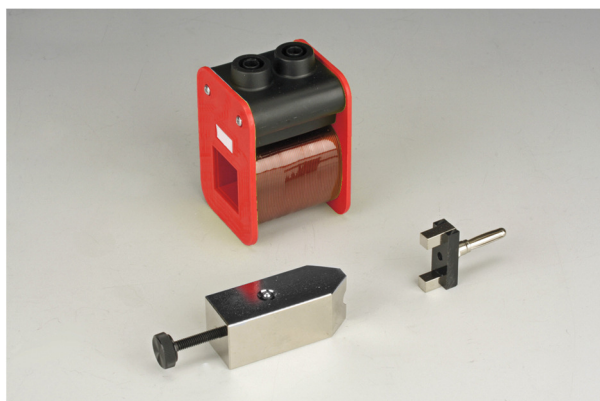
The launcher consists of

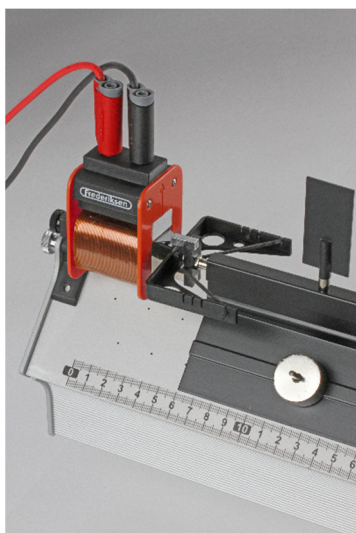
- Coil 1600 windings (462530)
- Iron core with finger screw
- Keeper with plug (195590)

Equipment needed

The iron core is designed to be fixed at the end of our air track (195050, 195000 or 195010). The keeper fits the carts of the air track.

The coil is placed on the iron core and is normally connected to a power supply through our switch box (198510 or 198500).

If an initial velocity is required, you can place one of the rubber band forks from the air track as shown.

Power supply requirements are 12 V and 0.5 A which leaves us with a lot of options (364000, 363000, 361600, 361800 ...).

In combination with the switch box, the electric launcher allows for reproducible releases of the air track cart – while simultaneously sending an electric start pulse to an electronic timer (200250).

Power supply

Normally, use a voltage of 12 V. The “8 V” text on the switch box does not apply for this version of the launcher.

Warning

Only turn on the current through the coil when it is actually required to hold the cart – i.e. shortly before launch.

However, the coil can without problems be connected to 12 V for up to an hour.

If a higher voltage than 12 V is used, the coil must be allowed to cool off between activations; at 24 V, keep a 40 seconds break for every 10 second the coil is turned on.

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© Frederiksen Scientific A/S