Beskrivelse

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2855.70 AC

Description

This spectral tube holder can be used with spectral tubes (Geissler tubes) between 224 and 299 mm long with a diameter of up to 15.5 mm.

The built-in power supply provides an ignition voltage of up to 6 kV. The operating voltage is significantly lower and is controlled by the variable current limiter. The current can be adjusted between 2 and 8 mA by means of a button on the back.

The tube is easily installed by inserting its lower end in the lower spring-loaded holder, press down and then put the top one in place.

Always turn off the power at the rear of the unit before replacing the tube, and let the tube cool before handling it. Do not turn the power back on until the new pipe is installed in place.

Never attempt to touch the electrodes in the bottom and top!

Current adjustment

The knob at the back of the cabinet is provided with a scale to read off the current in milliamps.

Spectral tubes have a limited life time. The life time is shortened by high operating temperatures, and the temperature increases with the current through the tube. Letting the tube cool down between measurements prolongs the lifetime significantly. For instance, you can alternate between 1 minute of operation and 5 minutes of cooling – in case it proves necessary to run the tube at the full 8 mA.

If the current is turned down too much, the tube may get unstable (blink) or eventually fail completely to turn on.

Technical data

L x W x H: 90 x 90 x 330 mm

Power consumption: 50 W

Mains voltage: 230 V, 50 Hz

Fuse: 315 mA, Slow

Accessories

The following spectral tubes are currently in stock:

2850.00 Ne	2550.10 Hg
2850.20 H ₂	2850.25 H ₂ Balmer
2850.30 He	2850.40 Ar
2850.50 O ₂ (IR)	2850.60 Kr
2850.70 N ₂	2851.10 H ₂ O
2851.20 CO ₂	2851.30 Xe
2851.50 I ₂	2851.80 Na and He

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© A/S Søren Frederiksen, Ølgod